

MiSeq Operating Software 1.0.40

Release Notes

For MiSeqDx™ Instruments

December 19, 2014

Introduction

These release notes describe all changes to MiSeq Operating Software for the MiSeqDx since the package containing MiSeq Operating Software 1.0.37.

This update is required for a diagnostic instrument. This update resolves an issue that could potentially lead to incorrect assignment of results to samples.

This update can be installed by the user. Follow the procedure in the customer install notes included in PCN0115.

The software package includes:

- MiSeq Operating Software 1.0.40

This release does not include changes to other software on the MiSeqDx instrument, including MiSeq Reporter, or to software run on another computer, such as Illumina Worklist Manager or Illumina User Management Software.

I. MiSeq Operating Software 1.0.40

NEW FEATURES:

- None

DEFECT REPAIRS:

- Fixed an issue with the handling of sample sheet file names. Previously, the inclusion of a period in the file name would result in the software attempting to load a sample sheet with a truncated file name. The software correctly handles periods in file names.
- Changed the default option for sending instrument health data to disabled.

KNOWN ISSUES:

- None